



**C E** Revised 12 Sept. 2011 rm (Vers. 8.1)

This kit is intended for Research Use Only.

Not for use in diagnostic procedures.

Please use only the valid version of the package insert provided with the kit

### INTENDED USE

Enzyme immunoassay for determination of melatonin sulfate (synonyms: 6-Hydroxymelatonin Sulfate, 6-Sulfatoxymelatonin) in human urine.

### TEST PRINCIPLE

Solid phase enzyme-linked immunosorbent assay (ELISA) based on the competition principle.

An unknown amount of antigen present in the sample and a fixed amount of enzyme labelled antigen compete for the binding sites of the antibodies coated onto the wells. After incubation the wells are washed to stop the competition reaction. After the substrate reaction the intensity of the developed color is inversely proportional to the amount of the antigen in the sample. Results of samples can be determined directly using the standard curve.

### WARNINGS AND PRECAUTIONS

- 1. For research use only. For professional use only.
- 2. Before starting the assay, read the instructions completely and carefully. Use the valid version of the package insert provided with the kit. Be sure that everything is understood.
- 3. In case of severe damage of the kit package please contact DRG or your supplier in written form, latest one week after receiving the kit. Don't use damaged components in test runs, but keep safe for complaint related issues.
- 4. Obey lot number and expiry date. Don't mix reagents of different lots. Don't use expired reagents.
- 5. Follow good laboratory practice and safety guidelines. Wear lab coats, disposable latex gloves and protective glasses where necessary.
- 6. Reagents of this kit containing hazardous material may cause eye and skin irritations. See MATERIALS SUPPLIED and labels for details. Material Safety Data Sheets for this product are available upon request.
- 7. Chemicals and prepared or used reagents have to be treated as hazardous waste according to national biohazard and safety guidelines or regulations.
- 8. Avoid contact with Stop solution. It may cause skin irritations and burns.

### STORAGE AND STABILITY

The kit is shipped at ambient temperature and should be stored at 2-8°C. Keep away from heat or direct sun light. The storage and stability of specimen and prepared reagents is stated in the corresponding chapters.

The microtiter plate is stable up to the expiry date of the kit in the broken, but tightly closed bag when stored at 2–8°C.





# DI CE Revised 12 Sept. 2011 rm (Vers. 8.1)

USA: RUO

### SPECIMEN COLLECTION AND STORAGE

### Urine

It is possible to use spontaneous as well as 24 h urine. The total volume of urine excreted during a 24 h period should be collected and mixed in a single bottle. Preservation is not necessary. Determine total volume for calculation of results. **Mix and centrifuge samples before use in the assay.** 

Storage:	2-8°C	≤ -20°C (Aliquots)	Keep away from heat or direct sun light. Avoid repeated freeze-thaw cycles.
Stability:	4 d	15 y	For more details see: Griefahn et al. (2001).

### MATERIALS SUPPLIED

Quantity	Symbol	Component		
1 x 12x8	MTP	<b>Microtiter Plate,</b> Break apart strips. Coated with anti-rabbit IgG (goat, polyclonal).		
1 x 5 mL	ANTISERUM	Melatonin Sulfate Antiserum; Ready to use.		
I X J IIIL	ANTISEKUM	Contains: Antiserum (rabbit), Tris buffer, 0.01 % Thimerosal		
	ENZCONJ CONC	Enzyme Conjugate Concentrate (40x)		
1 x 0.2 mL		Contains: Melatonin Sulfate, conjugated to peroxidase, phosphate buffer, 0.01 % Thimerosal.		
		Standard A-G; Ready to use.		
1 x 7 x 0.1 mL	CAL A-G	0; 1.7; 5.2; 15.6; 46.7; 140; 420 ng/mL		
		0; 5.2; 15.9; 47.6; 142; 427; 1281 nmol/L		
		Contains: Melatonin Sulfate, Tris buffer, 0.01 % Thimerosal.		
1 x 2 x	CONTROL 1+2	Control 1+2, Ready to use. Contains: 0.02 % Thimerosal.		
0.1 mL	CONTROL 1+2	Concentrations / acceptable ranges see QC certificate.		
1 x 60 mL	ASSAYBUF	Assay Buffer, Ready to use.		
T X 00 IIIL		Red colored. Contains: Tris buffer, BSA, 0.01 % Thimerosal.		
1 x 50 mL	WASHBUF	Wash Buffer Concentrate (20x)		
1 X JO IIIL	CONC	Contains: phosphate buffer, Tween, 0.1 % Thimerosal.		
1 x 12 mL	TMB SUBS	TMB Substrate Solution, Ready to use.		
		Contains TMB, buffer, stabilizers.		
1 x 12 mL	TMD STOP	TMB Stop Solution		
	TMB STOP	Ready to use. 1 M $H_2SO_4$ .		
3 x	FOIL	Adhesive Foil		





USA: RUO

# **C €** Revised 12 Sept. 2011 rm (Vers. 8.1)

### MATERIALS REQUIRED BUT NOT SUPPLIED

- 1. Micropipettes (Multipette Eppendorf or similar devices, < 3% CV). Volumes: 10; 50; 100; 1000 μL
- 2. Round-bottom polystyrene test tubes (12 x 75 mm)
- 3. Rack for test tubes
- 4. Orbital shaker (500 rpm)
- 5. Vortex mixer
- 6. 8-Channel Micropipettor with reagent reservoirs
- 7. Wash bottle, automated or semi-automated microtiter plate washing system
- 8. Microtiter plate reader capable of reading absorbance at 450 nm (reference wavelength 600-650 nm)
- 9. Bidistilled or deionised water
- 10. Paper towels, pipette tips and timer

### **PROCEDURE NOTES**

- 1. Any improper handling of samples or modification of the test procedure may influence the results. The indicated pipetting volumes, incubation times, temperatures and pretreatment steps have to be performed strictly according to the instructions. Use calibrated pipettes and devices only.
- 2. Once the test has been started, all steps should be completed without interruption. Make sure that required reagents, materials and devices are prepared ready at the appropriate time. Allow all reagents and specimens to reach room temperature (18-25 °C) and gently swirl each vial of liquid reagent and sample before use. Mix reagents without foaming.
- 3. Avoid contamination of reagents, pipettes and wells/tubes. Use new disposable plastic pipette tips for each component and specimen. Do not interchange caps. Always cap not used vials. Do not reuse wells/tubes or reagents.
- 4. Some components contain  $\leq 250 \ \mu$ L solution. Take care that the solution is completely on the bottom of the vial before opening.
- 5. It is advised to determine samples in duplicate to be able to identify potential pipetting errors.
- 6. Use a pipetting scheme to verify an appropriate plate layout.
- 7. Incubation time affects results. All wells should be handled in the same order and time sequences. It is recommended to use an 8-channel Micropipettor for pipetting of solutions in all wells.
- 8. Microplate washing is important. Improperly washed wells will give erroneous results. It is recommended to use a multichannel pipette or an automatic microplate washing system. Do not allow the wells to dry between incubations. Do not scratch coated wells during rinsing and aspiration. Rinse and fill all reagents with care. While rinsing, check that all wells are filled precisely with Wash Buffer, and that there are no residues in the wells.
- 9. Humidity affects the coated wells/tubes. Do not open the pouch until it reaches room temperature. Unused wells/tubes should be returned immediately to the resealed pouch including the desiccant.





# الکلات (( Revised 12 Sept. 2011 rm (Vers. 8.1)

### **PRE-TEST SETUP INSTRUCTIONS**

### Preparation of lyophilized or concentrated components

Dilute/ dissolve	Component		Diluent	Rela- tion	Remarks	Storage	Stability
15 mL	Wash Buffer	ad 300 mL	bidist. water	1:20	Resolve crystals at 18 – 25 °C.	2-8°C	4 w
50 µL	Enzyme Conjugate	with 2 mL	Assay Buffer	1:41	Prepare freshly and use only once.	18 – 25°C	30 min

### **Dilution of Standards, Controls and Urine Samples**

- Pipette 10 µL of each Standard, Control and urine sample into polystyrene, polypropylene or glass tubes. Avoid 1. direct sun light.
- Pipette 500 µL of Assay Buffer into each tube. Vortex. 2.

Samples containing concentrations higher than the highest standard have to be further diluted with Assay Buffer.

### **TEST PROCEDURE**

- 1. Pipette 50 µL of each diluted Standard, diluted Control and diluted sample into the respective wells of the Microtiter Plate.
- 2. Pipette **50** µL of freshly prepared **Enzyme Conjugate** into each well.
- 3. Pipette 50 µL of Melatonin Sulfate Antiserum into each well.
- Cover plate with adhesive foil. Incubate 2 h at RT (18-25°C) on an orbital shaker (500 rpm). 4.
- Remove adhesive foil. Discard incubation solution. Wash plate 4 x with 250  $\mu$ L of diluted Wash Buffer. Remove 5. excess solution by tapping the inverted plate on a paper towel.
- 6. For adding of Substrate and Stop Solution use, if available, an 8-channel Micropipettor. Pipetting should be carried out in the same time intervals for Substrate and Stop Solution. Use positive displacement and avoid formation of air bubbles.
- 7. Pipette 100 µL of TMB Substrate Solution into each well.
- 8. Incubate 30 min at RT (18-25°C) on an orbital shaker (500 rpm).
- Stop the substrate reaction by adding 100  $\mu$ L of TMB Stop Solution into each well. Briefly mix contents by gently 9. shaking the plate.
- 10. Measure optical density with a photometer at 450 nm (Reference-wavelength: 600-650 nm) within 60 min after pipetting of the Stop Solution.

### **QUALITY CONTROL**

The test results are only valid if the test has been performed following the instructions. Moreover the user must strictly adhere to the rules of GLP (Good Laboratory Practice) or other applicable standards/laws. All kit controls must be found

USA: RUO



# ىك ( ( Revised 12 Sept. 2011 rm (Vers. 8.1)

within the acceptable ranges as stated at the QC certificate. If the criteria are not met, the run is not valid and should be repeated. Each laboratory should use known samples as further controls. It is recommended to participate at appropriate quality assessment trials.

In case of any deviation the following technical issues should be proven: Expiration dates of (prepared) reagents, storage conditions, pipettes, devices, incubation conditions and washing methods.

### **CALCULATION OF RESULTS**

The obtained OD of the standards (y-axis, linear) are plotted against their concentration (x-axis, logarithmic) either on semi-logarithmic graph paper or using an automated method. A good fit is provided with cubic spline, 4 Parameter Logisitcs or Logit-Log.

For the calculation of the standard curve, apply each signal of the standards (one obvious outlier of duplicates might be omitted and the more plausible single value might be used).

The concentration of the samples can be read from the standard curve.

The initial dilution has been taken into consideration when reading the results from the graph. Results of samples of higher predilution have to be multiplied with the dilution factor.

Samples showing concentrations above the highest standard have to be diluted as described in PRE-TEST SETUP INSTRUCTIONS and reassayed.

Calculate the 24 h excretion for each urine sample:  $\mu g/24h = \mu g/L \times L/24h$ 

Conversion: Melatonin Sulfate  $(ng/mL) \times 3.05 = nmol/L$ 

### **Typical Calibration Curve**

(Example. Do not use for calculation!)

Standar d	Melatonin Sulfate (ng/mL)	Mean OD	OD/OD <sub>max</sub> (%)
А	0.0	1.805	100.0
В	1.7	1.741	96.5
С	5.2	1.536	85.1
D	15.6	1.185	65.7
Е	46.7	0.773	42.8
F	140	0.341	18.9
G	420	0.164	9.1

### **PRODUCT LITERATURE REFERENCES**

1. Kunz et al. Melatonin in s with Reduced REM Sleep Duration: Two Randomized Controlled Trials. Journal of Clinical Endocrinology & Metabolism, 89(1): 128-134 (2004)



USA: RUO







USA: RUO

### Revised 12 Sept. 2011 rm (Vers. 8.1)

- 1. Kondo, M. Tokura H., Wakamura T., Hyun KJ., Tamotsu S., Morita T., Oishi T., Influence of twilight on diurnal variation of core temperatuer, its nadir, and urinary 6-hydroxymelatonin sulfate during nocturnal sleep and morning drowsiness. Coll. Antropol. 33 (2009) 1: 193-199
- 2. Radwan P., Skrzydlo-Radomanska B., Radwan-Kwiatek K., Burak-Czapiuk B., Strzemecka J. Is melatonin involved in the irritable bowel syndrome? Journal of Physiology and Pharmacology (2009), Suppl 3, 67-70
- 3. Mahlberg R. et al. Normative data on the daily profile of urinary 6-sulfatoxymelatonin in healthy subjects between the ages of 20 and 84. Psychoneuroendocrinology (2006) 31, 634-641
- 4. Kunz et al. Melatonin in Patients with Reduced REM Sleep Duration: Two Randomized Controlled Trials. Journal of Clinical Endocrinology & Metabolism, 89(1): 128-134 (2004)
- 5. Mahlberg et al. Melatonin and age. Somnologie 7 (Suppl. 1): 9-88 (2003)
- 6. Griefahn et al. Melatonin production during childhood and adolescence: a longitudinal study on the excretion of urinary 6-hydroxymelatonin sulfate. J. Pineal Res. 34:26-31 (2003)
- 7. Bortkiewicz et al. The excretion of 6-hydroxymelatonin sulfate in healthy young men exposed to electromagnetic fields emitted by cellular phone – an experimental study. Neuroendocrinology Letters 23(suppl 1):88-91 (2002)
- 8. Kunz et al. Exogenous Melatonin in Periodic Limb Movement Disorder: An Open Trial and a Hypothesis. SLEEP, Vol. 24, No. 2, 183-187, (2001).
- 9. Griefahn et al. Long-Term Stability of 6-Hydroxymelatonin Sulfate in 24-h Urine Samples Stored at -20°C. Endocrine, vol. 15, no. 2, 199-202, July (2001).
- 10. Kunz D et al. Melatonin as a therapy in REM sleep behavior disorder patients: an open-labeled pilot study on the possible influence of melatonin on REM-sleep regulation. Movement Disorders, 14: 507-511 (1999)
- 11. Pfluger DH, Minder CE. Effects of exposure to 16.7 Hz magnetic fields on urinary 6-hydroxymelatonin sulfate excretion of Swiss railway workers. J. Pineal Res., 21: 91-100 (1996)

Version 2011-08-04~rm